

# **Technical Manual**

FluoroDazzle HRP Detection Reagent

**Catalogue Code: BA0173** 

Pack Size: 500 assay

**Research Use Only** 



## DESCRIPTION

The Assay Genie FluoroDazzle HRP Detection Reagent is a stable 10-Acetyl-3,7-dihydroxyphenoxazine (also called Ampliflu Red) liquid substrate reagent formulated for detecting horseradish peroxidase HRP activity, e.g. in ELISA. This reagent produces a highly fluorescent product ( $\lambda_{ex/em} = 530/590$ nm) when reacted with H<sub>2</sub>O<sub>2</sub> by a peroxidase. The fluorescent product can also be visualized under a light source ( $530 \pm 20$  nm).

### CONTENTS:

FluoroDazzle HRP Detection Reagent	25 mL Reagent
	100 $\mu L$ 3% Stabilized $H_2O_2$

The provided reagent is sufficient for 500 assays (96-well plate format).

## Instructions for use

Prior to assay, bring reagent to room temperature. Add 1  $\mu$ L 3% hydrogen peroxide per 1 mL HRP Reagent. For ELISA, use 50  $\mu$ L Reagent/H<sub>2</sub>O<sub>2</sub> per well in 96-well plate assays or 20  $\mu$ L per 384-well plate. Tap plate briefly, and incubate 30 min at room temperature. If desired, stop the reaction with 10 mM KOH (56 mg KOH/100 mL H<sub>2</sub>O). Read fluorescence at  $\lambda_{\text{ex/em}} = 530/590$ nm.

**Handling:** this product is shipped on ice. For long-term storage, keep reagent at -20°C. Shelf life of 1 year after receipt. **Bulk order** of any size (>125 mL) is available. Please contact info@assaygenie.com.

## LITERATURE

1. Towne V et al (2004). Complexities in horseradish peroxidase-catalyzed oxidation of dihydroxyphenoxazine derivatives: appropriate ranges for pH values and hydrogen peroxide concentrations in quantitative analysis. Anal Biochem. 334(2):290-6.

2. Li W, Jin W. (2006). Measurement of peroxidase activity in single neutrophils by combining catalyzed-enzyme reaction and epi-fluorescence microscopy. Talanta. 70(2):251-6.

#### PRECAUTIONS

Reagents are for research use only. Normal precautions for laboratory reagents should be exercised while using the reagents.

Avoid contact and inhalation. Standard laboratory safety procedures should be followed when handling this product. Safety procedures include wearing OSHA approved safety glasses, gloves and protective clothing.

If eye or skin contact occurs, immediately wash affected area with soap and copious amounts of water for at least 15 minutes. Assure adequate flushing by separating the eyelids with fingers. Call a physician to seek medical advice.

If swallowed, wash out mouth with water provided person is conscious and call a physician to seek medical advice.

If inhaled, move individual to fresh air and call a physician to seek medical advice.

#### TRANSPORT INFORMATION:

DOT (US):Not regulatedUN Number:Not regulatedClass:Not regulatedPackaging Group:Not regulated

Learn more at AssayGenie.com



**Contact Details** 

Dublin 2, Ireland

Email: info@assaygenie.com

Web: www.assaygenie.com

Technical Support: <u>Techsupportassaygenie.com</u>

Copyright © 2017 ReagentBio, All Rights Reserved. All information / detail is correct at time of going to print.

Learn more at AssayGenie.com