



# Murine IFN-gamma ELISpot Pair

**SKU:** MODC0179

Instructions for use

10x96 tests

**For research use only**

**Table of Contents**

1. Intended use ..... 2

2. Introduction ..... 2

2.1. Summary ..... 2

2.2. Principle of the method ..... 3

3. Reagents provided (Contents shown for 10x96 tests format) ..... 5

4. Materials/Reagents required but not provided ..... 5

5. Storage Instructions ..... 5

6. Safety & Precautions for use ..... 5

7. Reagent Preparation ..... 7

7.1. 1X Phosphate Buffered Saline (PBS) (Coating Buffer) ..... 7

7.2. Skimmed milk in PBS 1X solution (Blocking Buffer) ..... 7

7.3. 1% BSA PBS Solution (Dilution Buffer) ..... 7

7.4. 0.05% Tween PBS Solution (Wash Buffer) ..... 7

7.5. 35% Ethanol (PVDF Membrane Activation Buffer) ..... 7

7.6. Capture Antibody ..... 7

7.7. Detection Antibody ..... 7

7.8. Streptavidin conjugate ..... 8

7.9. Substrate buffer ..... 8

8. Sample and Control Preparation ..... 8

8.1. Cell Stimulation ..... 8

8.2. Positive Assay Control, mIFN-gamma production ..... 8

8.3. Negative Assay Control ..... 8

8.4. Sample ..... 8

9. Method ..... 9

10. Performance Characteristics ..... 10

10.1. Reproducibility and Linearity ..... 10

# Murine IFN-gamma ELISpot Pair

## 1. Intended use

Assay Genie **ELISpot** is a highly specific immunoassay for the analysis of cytokine and other soluble molecule production and secretion from T-cells at a single cell level in conditions closely comparable to the *in-vivo* environment with minimal cell manipulation. This technique is designed to determine the frequency of cytokine producing cells under a given stimulation and the comparison of such frequency against a specific treatment or pathological state. The ELISpot assay constitutes an ideal tool in the investigation of Th1 / Th2 responses, vaccine development, viral infection monitoring and treatment, cancerology, infectious disease, autoimmune diseases and transplantation.

Utilising sandwich immuno-enzyme technology, Assay Genie ELISpot assays can detect both secreted cytokines and single cells that simultaneously produce multiple cytokines. Cell secreted cytokines or soluble molecules are captured by coated antibodies avoiding diffusion in supernatant, protease degradation or binding on soluble membrane receptors. After cell removal, the captured cytokines are revealed by tracer antibodies and appropriate conjugates.

**This kit has been configured for research use only and is not to be used in diagnostic procedures.**

## 2. Introduction

### 2.1. Summary

Different populations of T-cells secrete differing patterns of cytokines that ultimately lead to different immune responses. IFN-gamma production is a key function of Th1, CD8<sup>+</sup> CTLs and also NK cells. IFN-gamma is a cytokine critical for cell mediated immunity against viral and intracellular bacterial infections and is involved in the inflammatory response following secretion via macrophage activation and stimulation of antibody secretion. IFN-gamma is the hallmark effector cytokine of Th1 and therefore is an excellent marker for identifying a host response to intracellular pathogens.

IFN-gamma is produced during infection by T cells of the cytotoxic/suppressor phenotype (CD8) and by a subtype of helper T cells, the Th1 cells. Th1 cells secrete IL-2, IL-3, TNF $\alpha$  and IFN-gamma, whereas Th2 cells mainly produce IL-3, IL-4, IL-5, and IL-10, but little or no IFN-gamma (1). IFN-gamma preferentially inhibits the proliferation of Th2 but not Th1 cells, indicating that the presence of IFN-gamma during an immune response will result in the preferential proliferation of Th1 cells (2).


In addition, IFN-gamma has several properties related to immunoregulation. IFN-gamma is a potent activator of mononuclear phagocytes (3), and activates macrophages to kill tumor cells by releasing reactive oxygen intermediates and TNF $\alpha$  (4). IFN-gamma induces or augments the expression of MHC antigens on macrophages, T and B cells and some tumor cell lines (5). On T and B cells IFN-gamma promotes differentiation. It enhances proliferation of activated B cells and can act synergistically with IL-2 to increase immunoglobulin light-chain synthesis (6, 7).

The role of IFN-gamma as a disease marker has been demonstrated for a number of different pathological situations including, viral infection (8), Autoimmune disease (9), transplant rejection (10), Diabetes (5) and allergy (11).


## 2.2. Principle of the method

A capture antibody highly specific for the analyte of interest is coated to the wells of a PVDF bottomed 96 well microtiter plate either during kit manufacture or in the laboratory. The plate is then blocked to minimise any non-antibody dependent unspecific binding and washed. Cell suspension and stimulant are added and the plate incubated allowing the specific antibodies to bind any analytes produced. Cells are then removed by washing prior to the addition of Biotinylated detection antibodies which bind to the previously captured analyte. Enzyme conjugated streptavidin is then added binding to the detection antibodies. Following incubation and washing substrate is then applied to the wells resulting in coloured spots which can be quantified using appropriate analysis software or manually using a microscope.


**1. 96-PVDF bottomed-well plates are first treated with 35% ethanol and then coated with capture antibody.**

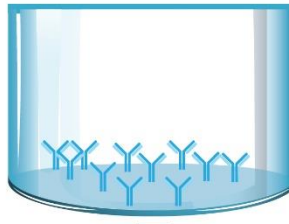


**2. Cells are incubated in the presence of the stimulating agent. Upon stimulation they release cytokines which bind to the capture antibodies.**

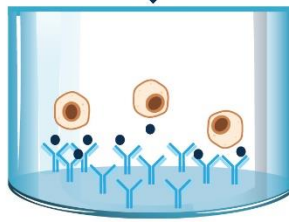

**3. Cells are removed by washing. Detection antibody is added.**

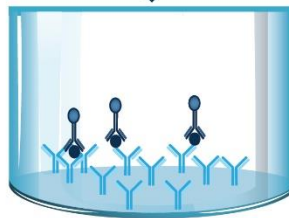


**4. Any excess unbound detection antibodies are removed by washing. Streptavidin-alkaline phosphatase is added.**

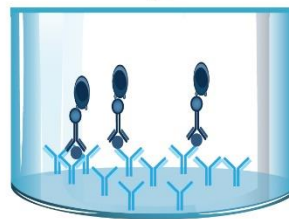





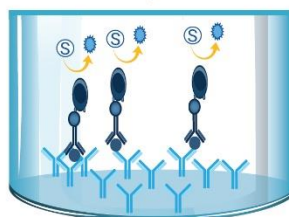

**5. Any excess unbound Strep-AP is removed by washing. BCIP/NBT is added.**

**BCIP/NBT is reduced by alkaline phosphatase to produce a precipitate observed as blue/purple spots.**

**One spot corresponds to one single producing cell.**



-  Capture antibody
-  Cytokines
-  Cell
-  Biotinylated detection antibody
-  Streptavidin-alkaline phosphatase conjugated
-  Substrate product

### 3. Reagents provided (Contents shown for 10x96 tests format)

- Capture Antibody (2 vials of 0.5 ml). The antibody is supplied sterile and does not contain preservative. We strongly advise sterile pipetting.
- Biotinylated detection antibody (2 vials, lyophilised)

### 4. Materials/Reagents required but not provided

- Miscellaneous laboratory plastic and/or glass, if possible sterile
- Streptavidin-conjugated (e.g. Streptavidin-alkaline phosphatase)
- Bovine Serum Albumin (BSA)
- Substrate solution (e.g. BCIP/NBT)
- Ethanol
- Cell culture reagents (e.g. RPMI-1640, L-glutamine, FCS)
- Cell stimulation reagents (PMA, Ionomycin)
- CO<sub>2</sub> incubator
- Tween 20
- Phosphate Buffered Saline (PBS)
- 96 well PVDF bottomed plates (we recommended Millipore plates catalogue # MSIPN4510, MSIPS4510 and M8IPS4510)

### 5. Storage Instructions

Store reagents between 2 and 8°C. Immediately after use remaining reagents should be returned to cold storage (2 to 8°C). Expiry of the components is stated on box front label and can only be guaranteed if the components are stored properly, and if in the case of repeated use of one component, the reagent is not contaminated by the first handling.

### 6. Safety & Precautions for use

- For **research use only** not to be used as a diagnostic test.
- Handling of reagents, blood specimens, PBMC, human cell lines should be in accordance with local safety procedures, e.g. CDC/NIH Health manual : "Biosafety in Microbiological and Biomedical Laboratories" 1984.
- Do not eat, drink, smoke or apply cosmetics where kit reagents are used.
- Do not pipette by mouth.
- When not in use, kit components should be stored refrigerated or frozen as indicated on vials or bottles labels.
- All reagents should be warmed to room temperature before use.
- Cover or cap all reagents when not in use.
- Do not mix or interchange reagents between different lots.
- Do not use reagents beyond the expiration date of the kit.
- Use a clean disposable plastic pipette tip for each reagent, standard, or specimen addition in order to avoid cross contamination.
- Use a clean plastic container to prepare the washing solution.
- Thoroughly mix the reagents and samples before use by agitation or swirling.

- All residual washing liquid must be drained from the wells by efficient aspiration or by decantation followed by tapping the plate forcefully on absorbent paper. Never insert absorbent paper directly into the wells.
- When pipetting reagents, maintain a consistent order of addition from well-to-well. This will ensure equal incubation times for all wells.
- **Substrate buffer** is potentially carcinogenic and should be disposed of appropriately, caution should be taken when handling this reagent, always wear gloves. Follow the supplier's instructions.
- Follow incubation times described in the assay procedure.

## 7. Reagent Preparation

### 7.1. 1X Phosphate Buffered Saline (PBS) (Coating Buffer)

For 1 litre of 10X PBS weigh-out:

80g NaCl
2g KH <sub>2</sub> PO <sub>4</sub>
14.4g Na <sub>2</sub> HPO <sub>4</sub> ; 2H <sub>2</sub> O.

Add distilled water to 1 litre.

**Dilute the solution to 1X before use.**

Check the pH of the 1X solution and adjust to required pH : 7.4 +/- 0.1.

### 7.2. Skimmed milk in PBS 1X solution (Blocking Buffer)

For one non-sterile plate, dissolve 0.2 g of dry skimmed milk in 10 ml of PBS 1X. For one sterile plate, dilute 5 ml of liquid milk in 5 ml of PBS 1X.

### 7.3. 1% BSA PBS Solution (Dilution Buffer)

For one plate, dissolve 0.2 g of BSA in 20 ml of PBS 1X.

### 7.4. 0.05% Tween PBS Solution (Wash Buffer)

For one plate, dilute 50 µl of Tween 20 in 100 ml of PBS 1X.

### 7.5. 35% Ethanol (PVDF Membrane Activation Buffer)

For one plate, mix 3.5 ml of ethanol with 6.5 ml of distilled water.

### 7.6. Capture Antibody

*This reagent is supplied sterile, once opened keep the vial sterile or aliquot and store at -20°C. For optimal performance prepare the Capture Antibody dilution immediately before use.*

For one plate, dilute 100 µl of capture antibody in 10 ml of PBS 1X and mix well.

### 7.7. Detection Antibody

Reconstitute the lyophilised antibody with 0.55 ml of distilled water. Gently mix the solution and wait until all the lyophilised material is back into solution.

*If not used within a short period of time, reconstituted Detection Antibody should be aliquoted and stored at -20°C. In these conditions the reagent is stable for at least one year. For optimal performance prepare the reconstituted antibody dilution immediately prior to use.*

For one plate, dilute 100 µl of antibody into 10 ml of Dilution Buffer and mix well.



### 7.8. Streptavidin conjugate

Dilute in Dilution buffer according to the instructions of the supplier.

### 7.9. Substrate buffer

Use according to the instructions of the supplier.

## 8. Sample and Control Preparation

### 8.1. Cell Stimulation

Cells can either be stimulated directly in the antibody coated wells (Direct) or, first stimulated in 24 well plates or flask, harvested, and then plated into the coated wells (Indirect).

The method used is dependent on 1) the type of cell assayed 2) the expected cell frequency. When a low number of cytokine producing cells are expected it is also advised to test them with the direct method, however, when this number is particularly high it is better to use the indirect ELISpot method.

All the method steps following stimulation of the cells are the same whatever the method (direct/indirect) chosen.

### 8.2. Positive Assay Control, mIFN-gamma production

We recommend using the following polyclonal activation as a positive control in your assay.

Dilute mouse splenocytes in culture medium (e.g. RPMI 1640 supplemented with 2mM L-glutamine and 10% heat inactivated fetal calf serum) containing 0.5 µg/ml Concanavalin A and 2 ng/ml mL-2. Incubate 2 days (Indirect stimulation). Take off non adherent materials and harvest adherent cells with a cells scraper. Wash cells once.

Dilute cells in culture medium supplemented with 1 ng/ml PMA and 500 ng/ml ionomycin (Sigma, Saint Louis, MO). Distribute  $1 \times 10^5$  to  $2.5 \times 10^5$  cells per 100 µl in required wells of an antibody coated 96-well PVDF plates and incubate for 15-20 hours in an incubator.

For other stimulators incubation times may vary, depending on the frequency of cytokine producing cells, and should be optimised in each situation.

### 8.3. Negative Assay Control

Dilute mouse splenocytes in culture medium to give an appropriate cell number (same number of unstimulated cells as stimulated sample cells) per 100 µl with no stimulation.

### 8.4. Sample

Dilute mouse splenocytes in culture medium and stimulator of interest (i.e. Sample, Vaccine, Peptide pool or infected cells) to give an appropriate cell number per 100 µl.

Optimal assay performances are observed between  $1 \times 10^5$  and  $2.5 \times 10^5$  cells per 100  $\mu\text{l}$ .

Stimulators and incubation times can be varied depending on the frequency of cytokine producing cells and therefore should be optimised by the testing laboratory.

## 9. Method

Prepare all reagents as shown in section 7 and 8.

**Note: For optimal performance prepare the Streptavidin-AP dilution immediately prior to use.**

Assay Step		Details
1.	Addition	For PVDF membrane activation, add 25 $\mu\text{l}$ of 35% ethanol to every well
2.	Incubation	Incubate plate at room temperature (RT) for 30 seconds
3.	Wash	Empty the wells by flicking the plate over a sink & gently tapping on absorbent paper. Thoroughly wash the plate 3x with 100 $\mu\text{l}$ of PBS 1X per well
4.	Addition	Add 100 $\mu\text{l}$ of diluted <b>capture antibody</b> to every well
5.	Incubation	Cover the plate and incubate at 4°C overnight
6.	Wash	Empty the wells as previous and wash the plate once with 100 $\mu\text{l}$ of PBS 1X per well
7.	Addition	Add 100 $\mu\text{l}$ of <b>blocking buffer</b> to every well
8.	Incubation	Cover the plate and incubate at RT for 2 hours
9.	Wash	Empty the wells as previous and thoroughly wash 3x with 100 $\mu\text{l}$ of PBS 1X per well
10.	Addition	Add 100 $\mu\text{l}$ of <b>sample, positive and negative controls</b> cell suspension to appropriate wells providing the required concentration of cells and stimulant (cells may have been previously stimulated see section 8.)
11.	Incubation	Cover the plate and incubate at 37°C in a CO <sub>2</sub> incubator for an appropriate length of time (15-20 hours). <b>Note: do not agitate or move the plate during this incubation</b>
12.	Addition	Empty the wells and remove excess solution then add 100 $\mu\text{l}$ of Wash Buffer to every well
13.	Incubation	Incubate the plate at 4°C for 10 min
14.	Wash	Empty the wells as previous and wash the plate 3x with 100 $\mu\text{l}$ of Wash Buffer
15.	Addition	Add 100 $\mu\text{l}$ of diluted <b>detection antibody</b> to every well
16.	Incubation	Cover the plate and incubate at RT for 1 hour 30 min
17.	Wash	Empty the wells as previous and wash the plate 3x with 100 $\mu\text{l}$ of Wash Buffer
18.	Addition	Add 100 $\mu\text{l}$ of diluted <b>Streptavidin-conjugate</b> to every well
19.	Incubation	Cover the plate and incubate at RT following the supplier's instructions
20.	Wash	Empty the wells and wash the plate 3x with 100 $\mu\text{l}$ of Wash Buffer
21.	Wash	Peel of the plate bottom and wash both sides of the membrane 3x under running distilled water, once washing complete remove any excess solution by repeated tapping on absorbent paper.
22.	Addition	Add 100 $\mu\text{l}$ of ready-to-use <b>substrate buffer</b> to every well

23.	Development	Following the supplier's instructions, incubate the plate for <b>5-15 min</b> monitoring spot formation visually throughout the incubation period to assess sufficient colour development
24.	Wash	Empty the wells and rinse both sides of the membrane 3x under running distilled water. Completely remove any excess solution by gentle repeated tapping on absorbent paper

**Read Spots:** allow the wells to dry and then read results. The frequency of the resulting coloured spots corresponding to the cytokine producing cells can be determined using an appropriate ELISpot reader and analysis software or manually using a microscope.

*Note: spots may become sharper after overnight incubation at 4°C in the dark*

Plate should be stored at RT away from direct light, but please note that colour may fade over prolonged periods so read results within 24 hours.

## 10. Performance Characteristics

### 10.1. Reproducibility and Linearity

Intra-assay reproducibility and linearity were evaluated by measuring the spot development following the stimulation (PMA / Ionomycin) of 5 different mouse splenocytes cell concentrations, 12 repetitions. The data show the mean spot number, range and CV for the five cell concentrations.

Cells / well	n	Mean number of spots per well	Min	Max	CV%
100000	12	462	406	495	5.5%
50000 recommended	12	494	457	518	3.7%
25000 recommended	12	391	355	411	4.8%
12500	12	244	202	267	7.8%
6250	12	143	120	190	12.5%



**Contact Us:**

Assay Genie  
The Steelworks  
Dublin  
Ireland  
D01 KA00

**Email:** [info@assaygenie.com](mailto:info@assaygenie.com)

**Phone:** 00-353-1-887-9802